



## Rules for operating and cleaning the cryostat

2020-03-19 version 1

All users must attend an introduction meeting.

Shared by all

- Bottle with detergent & brush for cleaning the sample stubs
- Bottle with distilled water

Company owned

- Sectioning equipment (knives, cryomold etc)
- EtOH for cleaning
- Brushes

Cleaning after use

- Remove the knife and clean the chamber from sectioning residue
- Wipe clean with 96% EtOH in the chamber
- Clean the brushes with 96% EtOH
- Clean the sample stubs by scrubbing them with water and detergent, rinse with distilled water.
- Discard waste into waste box

SmiLe will change the anti-roll glass plate 1-2 times / year.

Companies are welcome to bring their own plate.

SmiLe provides waste box and disposal thereof.