
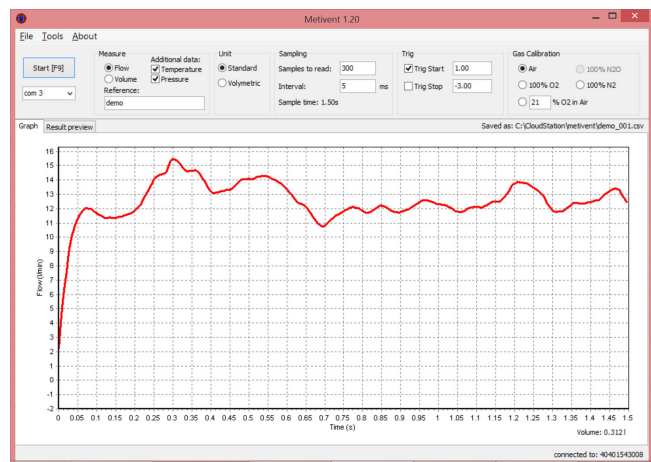


Metivent Software

A Companion to TSI Flowmeters

Metivent is a Windows application for recording air flow measurements from TSI flowmeters (e.g. TSI4040). A standard Windows computer with the installed Metivent software connects to the flowmeter via a USB-serial converter. The acquisition of data from the flowmeter is then controlled by the software. Once started, the graph is updated continuously during measurement.

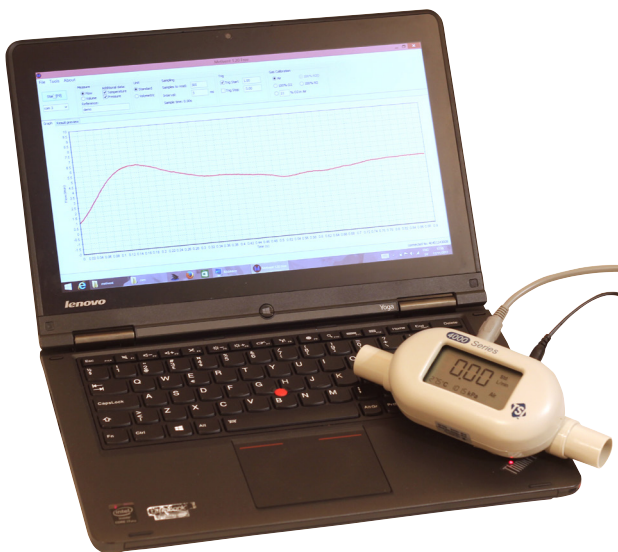

The software can be used in several modes and configurations.

- Start trig (wait to start until the flow has reached the start trigger level)
- End trig (sample data until stop trigger condition)
- Gas calibration
- Number of samples to collect if not using trig
- Standard or volumetric flow
- Sampling rate 1-1000/s

Measurements can be conveniently saved to CSV files and opened in Excel for further data analyses and presentations. The instrument is pivotal when recording inhalations profiles for lung simulation studies (such as FIA's F-SIG 6300).

The file presents flow, pressure and temperature along with instrument data, e.g. serial number, calibration date and currently used settings.

	A	B	C	D	E
1	Sample Reference	20170105-sample1			
2					
3	Program Version	1.2			
4					
5	Instrument:				
6	Model	4040			
7	Serial/ID#	4040213904			
8	Calibration Date	2014-10-15			
9	Firmware	3.5			
10					
11	Settings:				
12	Measurement	Flow			
13	Read Temp	Yes			
14	Read Pressure	Yes			
15	Flow Unit	Standard			
16	Samples	300			
17	Sample interval (ms)	3			
18	Trig Start	NA			
19	Trig Stop	NA			
20	Gas Calibration	Air			
21					
22	Time(ms)	Flow (l/min)	Temp (C)	Pressure (kPa)	
23	0	0.01	27.01	102.22	
24	3	0	27.01	102.19	
25	6	0	27.01	102.29	
26	9	0	27.01	102.14	
27	12	0	27.01	102.09	
28	15	0	26.99	102.15	
29	18	0	26.98	102.25	
30	21	0	26.98	102.07	
31	24	0	26.98	102.22	
32	27	0	26.98	102.32	
33	30	0.01	26.98	102.14	
34	33	0.03	26.98	102.1	
35	36	0.03	26.98	102.29	

For contact and more information

Kjell Fransson
AB FIA
Managing Director
kjell.fransson@fia.se
Mobile: +46 701 45 54 83

Member of
MVIC
MEDICIN VALLEY
INHALATION CONSORTIUM
www.mvic.se



www.fia.se

AB FIA
An Engineering Company
with a Chemistry Profile

AB FIA Vinkelhaken 1 D
SE-247 32 Södra Sandby
Sweden • i@fia.se